



Bi-Test™ CD3 FITC - CD62L PE

Product: Anti-human T cell monoclonal I antibody and anti-human L-Selectin monoclonal antibody.

Description: The CD3 epitope is expressed on the epsilon chain of the CD3/T cell antigen receptor (TcR) complex. CD3 is expressed on 65-85% of thymocytes and has a mitogenic effect on peripheral blood T cells. Identification of human T cells expressing the 22-28,000 M. W. surface antigen. CD62L is the equivalent of the mouse MEL-14 with a molecular weight of 80 KDa. CD62L is on 60-75% of peripheral blood lymphocytes, Approximately 70-80% of T cells, 10% of thymocytes and most B cells. CD62L reacts with monocytes, granulocytes and a subset of NK cells. The selectins represent a family of adhesion molecules, expressed by leukocytes and endothelial cells that are involved in the regulation of leukocyte traffic. Human leukocyte L selectin is also called LAM-1, LECAM-1, or gp90MEL-14. This antibody to L selectin blocks greater than 95% of lymphocyte adhesion to peripheral lymph node high endothelial venules.

Isotype: Mouse IgG1 kappa (FITC) and Mouse IgG2a kappa (PE)

Clones: M2AB (CD3 FITC) and 1H3 (CD62L PE)

Applications: Monitoring of T cells subsets in peripheral blood; Characterization of subtypes of T cell leukemia's and lymphomas; Studies of AIDS/HIV virus infection; Analysis of CD3 complex related to the T cell antigen receptor; Study of Cell-adhesion molecules; Study of B cell functions.

Use: PBMC: Add 10 µl of MAB/10⁶ PBMC in 100 µl PBS. Mix gently and incubate for 15 minutes at 2^o to 8^o C. Wash twice with PBS and analyze or fix with 0.5% v/v of paraformaldehyde in PBS and analyze. WHOLE BLOOD: Add 10 µl of MAB/100 µl of whole blood. Mix gently and incubate for 15 minutes at room temperature 20^o C. Lyse the whole blood. Wash once with PBS and analyze or fix with 0.5% v/v of paraformaldehyde in PBS and analyze. See instrument manufacturer's instructions for Lysed Whole Blood and Immunofluorescence analysis with a flow cytometer or microscope.

Storage: Antibodies are supplied in PBS, 0.08% sodium azide and 0.2% protein carrier for FITC and PE. Antibodies should be stored at 4-8^oC. Monoclonal antibodies should not be frozen. Reagents are stable for the period shown on the vial label when stored properly.

Ordering Information:	Form	Vial Size	Catalog #
	Bi-Test™	50 Test	362Ls
	Bi-Test™	100 Test	362L

For research use only. Not for use in human diagnostics or therapeutics.

REFERENCES:

1. Knowles RW. Immunochemical analysis of the T cell-specific antigens. In : Reinhert EL, Haynes BF, Nadl LM and Bernstein ID. eds. Leukocyte Typing II, Human T Lymphocytes. New York, NY: Springer-Verlag; 1986:259.
2. Kurre R. Cluster Report:CD3. In:Knapp W, Dorken B, Gilks WR, Reiber EP, Schmidt RE, Stein H, and von dem Borne AEG Kr, eds. Leukocyte Typing IV, White cell Differentiation Antigens. Oxford, England: Oxford Press 1989:293.
3. Signal transduction via CD4,CD8 and CD28 in mature and immature thymocytes. Implications for thymic selection. Turka LA, Linsley PS, Paine R 3d, Schieven GI, Thompson GB, Ledbetter JA, J. Immunol. 1991 Mar :146(5): 1428-36.
4. T cell receptor/CD3-signaling induces death by apoptosis in human T cell receptor gamma delta + Tcells. Janssen O, Wesselborg S, Heckl-Ostreicher B, Pechhold K, Bender A, Schondelmaier S, Moldenhauer G, Kabelitz D I Immunol. 1991 Jan146(1):35-9.
5. Clonal analysis of human CD4-CD8-CD3- thymocytes highly purified from postnatal thymus Hori T, Spits H J. Immunol. 1991 Apr 146(7):2116-21.
6. Molecular cloning of the CD3 zeta subunit identifies a CD3 zeta-related product in thymus-derived cells Jin YJ, Claton LK, Howard FD, Koyasu S, Sieh M, Steinbrich R, Tarr GE, Reinherz EL. Proc Natl Acad Sci USA 1990 Ma: 87(9):3319-23.
7. Immunofluorescence Measurement in a Flow Cytometer using Low-Power Helium Neon Laser Excitation. Shapiro, H.M, Glazer, A.N., Christenson, L., Williams, J.M., and Strom, T. B. Cytometry 4,276, 1983.
8. Comparison of Helium Neon and Dye lasers for Excitation of Allophycocyanin. Loken, M.R., Kiej, J.F. and Kelly, K.,A. Cytometry 8, 96, 1987.
9. Identification of a specific glycoprotein ligand for P-selectin (CD62) on myeloid cells. J Cell Biol 1992 Jul;118(2):445-56. Moore KL; Stults NL; Diaz S; Smith DF; Cummings RD; Varki A; McEver RP
10. Cytoplasmic domain of P-selectin (CD62) contains the signal for sorting into the regulated secretory pathway. Mol Biol Cell 1992 Mar;3(3):309-21. Disdier M; Morrissey JH; Fugate RD; Bainton DF; McEver RP
11. Leukocytes roll on a selectin at physiologic flow rates: distinction from and prerequisite for adhesion through integrins. Cell 1991 May 31;65(5):859-73. Lawrence MB; Springer TA
12. Biochemical and immunohistochemical characteristics of CD62 and CD63 monoclonal antibodies. Expression of GMP-140 and LIMP-CD63 (CD63 antigen) in human lymphoid tissues. Virchows Arch B Cell Pathol 1991;61(4):269-77. Metzelaar MJ; Schuurman HJ; Heijnen HF; Sixma JJ; Nieuwenhuis HK
13. CD62/P-selection recognition of myeloid and tumor cell sulfatides. Cell 1991 Oct 4;67(1):35-44. Aruffo A; Kolanus W; Walz G; Fredman P; Seed B
14. Structure and chromosomal location of the gene for endothelial-leukocyte adhesion molecule 1. J Biol Chem 1991 Feb 5;266(4):2466-73. Collins T; Williams A; Johnston GI; Kim J; Eddy R; Shows T; Gimbrone MA Jr; Bevilacqua MP

For research use only. Not for use in human diagnostics or therapeutics.

Exalpa Biologicals, Inc., 86 Rosedale Rd. Watertown, MA 02472
Tel: 800.395.1137 or 617.924.3400, Fax: 866.924.5100 or 617.924.5100, Web:www.exalpa.com